RIDA[®] Sample Decolorant

Art. No.: R1699

In vitro Test

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Storage at 2 - 25 °C (35 - 77 °F)

R-Biopharm AG, Darmstadt, Germany

Tel.: +49 (0) 61 51 81 02-0 / Telefax: +49 (0) 61 51 81 02-20

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RIDA[®] Sample Decolorant

1. Intended use

RIDA[®] Sample Decolorant (patent pending EP 160 000 276.2) is used for the preparation of samples, such as:

- wine for RIDASCREEN[®] Histamin (enzymatic) (R1605)
- liver, feed for RIDASCREEN[®] Clenbuterol (R1711)

Thereby, RIDA[®] Sample Decolorant removes interfering substances by precipitation. Thereafter the samples can be analyzed according to the respective instructions for use or application note.

2. Reagents provided

The reagents of one package are sufficient to extract e.g. 200 wine samples, 400 liver samples or 200 feed samples.

Component	CAP colour	Format	Volume
Reagent 1	colorless	Ready-to-use	40 ml
Reagent 2	blue	Ready-to-use	40 ml
Reagent 3	green	Ready-to-use	20 ml

3. Warnings and precautions for users

Reagent 2 (Decolorant) is harmful to health. It should be worked under a chemical hood and skin contact should be avoided.

The solutions should only be used by trained laboratory employees.

This kit may contain further hazardous substances. For hazard notes on the contained substances please refer to the appropriate material safety data sheets (MSDS) for this product, available online at www.r-biopharm.com.

4. Storage instructions

Store at 2 - 25 °C / 35 - 77 °F.

Do not allow product to reach sewage system. Disposal of waste according to your national regulations.

No quality guarantee is accepted after the expiry date on the kit label.

5. Procedure

Bring reagents to room temperature before use. If crystals have formed dissolve them by shaking.

Perform sample preparation according to respective instructions for use or application note.

The data corresponds to our present state of technology and provides information on our products and their uses. R-Biopharm makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. Defective products will be replaced. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. R-Biopharm shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product

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are registered trademarks of R-Biopharm AG